

Maintenance Plan of Z-Arm Robotic Arm										
Position of Maintenance	Place of Maintenance	Once /Quarter	Once /Half a year	Once/ 1 year	Once/ 2years	Once/ 3 years	Methods of Maintenance	Time of Maintenance	Qty Per Set	Sale Price ( Unit Price )
Screw that is easily loosened by impact	Robot base fixing screw			√			Fastening	5minutes		
	Load mounting screw of robotic arm end	√					Fastening	5minutes		
	Routing of robotic arm end	√					Fastening	5minutes		
	Robot arm fixing screw		√				Fastening	10minutes		
	Rotary joint fixing screw of robotic arm end	√					Fastening	2minutes		
	Fixing screw of main body housing		√				Fastening	2minutes		
Is the cable loose or damaged?	Power socket of robotic arm		√				Fastening	20minutes		
	Communication cable of Robotic arm		√				Fastening	20minutes		
Whether the appearance is dirty, bumped, etc.; if so, please clean or replace it in time.	Surface of robotic arm			√			Wiping	30minutes		
	Power and control cable outer surface		√				Wiping	5minutes		
	Servo motor encoder battery replacement					√	Please contact the original factory for replacement if there is no power.		3	30
	Servo motor bearing					√	Please contact the original factory for replacement if there is damaged.		6	70

	Up and down motion towline				√		Please contact the original factory for replacement if there is damaged.		1	100
--	-------------------------------	--	--	--	---	--	---	--	---	-----

Huiling-tech Robotic Co., Ltd